

Primer & Probe reagent dilution mannul

Name	RT Primer	Scale	1nmol
Purification	PAGE	Attachment	1.5mL DNase/RNase-free ddH ₂ O
Format	Lyophilized		
Dilution	Add 100μL DNase/RNase-free ddH ₂ O to 10μM RT primer dilution		

Name	PCR primer set	Scale	2nmol
Purification	PAGE	Attachment	1.5mL DNase/RNase-free ddH ₂ O
Format	Lyophilized		
Dilution	Add 200μL DNase/RNase-free ddH ₂ O to 10μM primer set dilution		

Name	TaqMan probe	Scale	1nmol
Purification	HPLC	Attachment	1.5mL DNase/RNase-free ddH ₂ O
fluorescence Lable	5' FAM 3'BHQ1		
Format	Lyophilized and keep in dark		
Dilution	Add 100μL DNase/RNase-free ddH ₂ O to 10μM probe dilution		

Storage: Lyophilized powder could be stored at -20°C for 2 years, dilution at -20°C for 6 months to 1 year and at 4°C for 1 month.

Note: Please centrifuge the tube for 1 minute at 10,000–12,000×g before open it carefully. Add approximate DNase/RNase-free ddH₂O to the bottom of tube, then close the tube and vortex.

Contact us

Phone: 021-51320195-8002(Shanghai) 0512-86668828-8042(Suzhou)

E-mail: support@genepharma.com(Shanghai)

szsupport@genepharma.com(Suzhou)

<http://www.genepharma.com>



引物探针试剂稀释说明

类型	逆转录引物	规格	1nmol
纯化方式	PAGE	附送	1.5mL DEPC处理水
形态	减压离心干燥品，制品干燥于管底，为干粉状		
溶液配制	为配制成 10μM 溶液，需要在每一管中加入 100μL DEPC处理水		

类型	PCR 引物对	规格	2nmol
纯化方式	PAGE	附送	1.5mL DEPC处理水
形态	减压离心干燥品，制品干燥于管底，为干粉状		
溶液配制	为配制成 10μM 溶液，需要在每一管中加入 200μL DEPC处理水		

类型	TaqMan 探针	规格	1nmol
纯化方式	HPLC	附送	1.5mL DEPC处理水
荧光标记	5' 标记 FAM 3'标记 BHQ1		
形态	减压离心干燥品，制品干燥于管底，为干粉状，需避光保存		
溶液配制	为配制成 10μM 溶液，需要在每一管中加入 100μL DEPC处理水		

贮存条件：减压离心干燥品可于-20℃保存 2 年，溶液可于-20℃ 保存 6 个月到 1 年，可于4℃保存 1 个月；

注意：由于DNA oligo 呈很轻的干膜附在管壁上，打开时极易散失，所以开盖前应先离心10,000–12,000 $\times$ g 1 分钟，然后慢慢打开管盖，溶解时请加适量的DEPC 处理水后盖上管盖，振荡溶解。

联系我们

电话：021-51320195-8002(上海) 0512-86668828-8042(苏州)

E-mail: support@genepharma.com(上海)

szsupport@genepharma.com(苏州)

http://www.genepharma.com



B024-V005-20191028